

Whole Mount Retina IHC

Day 1

1. Fix eyes in **4% PFA** for 45-60"
2. Wash 3 x 10" in PBS
3. Dissect retinas
4. Block with **1% BSA/0.5% Tween 20** in PBS for 30" at rt
5. Wash 2x in **PBlec** (0.1 nM CaCl₂, 0.1 MgCl₂ mM, 0.1 MnCl₂ mM, and 1% Triton X-100 in PBS), incubate in last wash at least 30'.
6.
 - a. Incubate 500 ml o/n at 4°C with **Isolectin B4-biotin (1:10)** (SIGMA, L-2140) in PBlec, shaking.
 - b. Incubate 500 ml o/n at 4°C with Lectim-FITIC 1:50 (Sigma 107K4118) in PBlec, shaking. 2 minutes maximum velocity RT.

Day 2

1. Wash 5 x 10" in PBS at rt
2. Incubate o/n at 4°C with **Streptavidin-Alexa 488 (1:100)** centrifuge before use (Molecular Probes, S11223) and DAPI (1:1000) (Invitrogen, D3571) in 0.5% BSA/0.25% Tween 20 in PBS, shaking

Day 3

1. Wash 5 x 10" in PBS at rt
2. Mount in **ProLong Gold anti-fade** (Invitrogen, P36934)

**HAMSTER ANTI MOUSE CD31
MONOCLONAL ANTIBODY**

CATALOG NUMBER: MAB1398Z

LOT NUMBER: xxxxxx

QUANTITY: 100 µg

CONCENTRATION: 1 mg/mL

SPECIFICITY: React s with the 130-140 kDa protein, PECAM-1 found on mouse monocytes, platelets, granulocytes and endothelial cells. 2H8 effectively inhibits transmigration of activated polymophonuclear cells and monocytes across the endothelium. In a mouse model for acute peritonis - 2H8 blocks acute inflammation.

ISOTYPE: IgG

CLONE NAME : 2H8

APPLICATIONS: Flow Cytometry: 0.1 µg/mL
Blocking
Optimal working dilutions must be determined by end user.

FORMAT: Purified immunoglobulin - Ig fraction

PRESENTATION: Liquid in phosphate buffered saline containing no preservatives

STORAGE/HANDLING: Store at -20°C only. Avoid repeated freeze-thaw cycles.

REFERENCES:

1. Yong, X et al. (1993). *Proc. Natl. Acad. Sci.* **90**: 5569-5573
2. Bogen, S. et al. (1994). *J. Exp. Med.* **179**: 1059-1064.
3. Newman, P.J. et al. (1992). *J. Cell. Biol.* **119**: 239-246.

For research use only; not for use as a diagnostic.

□ 2003-2005: CHEMICON International, Inc. - By CHEMICON International, Inc. All rights reserved. No part of these works may be reproduced in any form without permissions in writing.

USA & Canada • Phone: +1(800) 437-7500 • Fax: +1 (951) 676-9209 • Europe +44
(0) 23 8026 2233
Australia +61 3 9839 2000 • Germany
+49-6192-207300 • ISO
Registered worldwide
www.chemicon.com •
custserv@chemicon.com •
techserv@chemicon.com

061704/mab1398z/ef/02/03/05/ft